Little Dipper[®] Processor

USER MANUAL

FOR RESEARCH USE ONLY

Warranty

SciGene warrants that the electrical instrument described in this manual shall be free of defects in materials and workmanship for a period of 12 months from date of delivery. This warranty does not cover removable baths or accessories. In the event of a defect during the warranty period, SciGene's limit of liability will be to provide replacement parts at no charge or, at its sole discretion, replace the product. The foregoing warranty is void in the event the unit was abused or modified or used in a manner inconsistent with its intended purpose. SciGene makes no other warranty, expressed or implied including warranties of merchantability and fitness for a particular purpose. In no event shall SciGene be liable for any direct, indirect, special, incidental or consequential damages or for any damages resulting from loss arising out of or in connection with the sale, use or performance of the product.

Copyright

Copyright ©2005-2023 SciGene Corporation. All rights reserved. SciGene and Little Dipper are trademarks of SciGene Corporation, Sunnyvale, CA. All other trademarks used in this manual are the property of their respective owners.

Table of Contents

I. SAFETY NOTICES	3
A. Intended Use	3
B. Instrument Safety.....	3
C. Symbols and Conventions	3
D. Warnings	4
E. Cautions.....	5
F. Spills and Reagent Disposal	6
G. Compliance.....	7
II. UNPACKING AND SET UP	8
A. Unpacking the Instrument	8
B. Parts Provided	8
C. Environmental Requirements	8
D. Installation and Location	8
E. Installing Baths	9
F. Installing Centrifuge Bucket Assemblies	9
III. USING YOUR LITTLE DIPPER PROCESSOR	10
A. Components, Controls and Accessories	10
B. Touch Screen and Power Connections.....	11
C. Using the Temperature Controller	12
D. Calibrating the Temperature Controller	12
IV. SYSTEM OPERATION	13
A. Using the Touchscreen	13
B. Using the Pause Button	13
C. Using the Emergency Stop (EMO)	13
D. Using a Bath Cover for Pre-Heating	14
E. Loading a Slide Rack	15
F. Running a Programmed Protocol	16
V. USING CUSTOM PROTOCOLS	17
A. Creating or Editing a Protocol	17
B. Deleting a Protocol.....	18
VI. MAINTAINING THE LITTLE DIPPER PROCESSOR	19
A. Powering Off	19
B. Cleaning Before Every Use	19
C. Cleaning After Every Use	19
D. Periodic Cleaning and Calibration	19
VII. TROUBLESHOOTING	20
VIII. SPECIFICATIONS	20
IX. ORDERING INFORMATION	20
X. DECLARATION OF CONFORMITY	21

I. SAFETY NOTICES

A. Intended Use

The Little Dipper Processor automates slide washing and drying for FISH (fluorescence in situ hybridization), chromosome and microarray analyses.

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B. Instrument Safety

Before operating the instrument, read the information in this section concerning hazards and potential hazards. Ensure that anyone involved with the instrument's operation is instructed in both general safety practices for laboratories and specific safety practices for the instrument.

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C. Symbols and Conventions

The following list is an illustrated glossary of the electrical symbols that are used on the **Little Dipper Processor**. Whenever such symbols appear on instruments, please observe appropriate safety measures.

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1. Electrical Symbols

This symbol indicates that this is a protected **ground terminal** that must be connected to earth ground before any other electrical connections are made to the instrument.

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CAUTION: This symbol alerts you to an **electrical hazard**. Please consult this User Manual for further information and proceed with caution.

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This symbol indicates the OFF position of the **main POWER** switch.

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This symbol indicates the ON position of the **main POWER** switch.

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D. Warnings

Failure to comply with the following warnings that are affixed to the product can lead to possible personal injury or death.

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E. Cautions

Failure to comply with the following cautionary statement affixed to the product may lead to possible personal injury.

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This symbol indicates the potential for a **pinch point**. Do not place fingers in the slot in the Z-axis tower.

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This symbol indicates the potential presence of a **hot surface**. Turn off the temperature controller and leave sufficient time for cooling or use heat protection prior to handling a bath or cleaning in this area.

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Do not touch the bottom **metal surface** of the bath during operation.

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Biological hazards can be associated with the materials used on this instrument. Potentially infectious waste and other hazardous waste must be treated in accordance with good laboratory practice guidelines.

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This symbol on the rear of the instrument indicates the presence of the fuse box. **Warning: For Continued Protection Against Fire, Replace Only with Same Type Rating of Fuse.** Always disconnect the power cord before attempting to replace the fuse.

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Always insert a **bath liner** and fill with buffer before turning on power to each bath.

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The instrument weighs approximately 59 lbs (26.5 kg). Use caution when **lifting** the unit to avoid personal injury.

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Do not under any circumstances lift the unit by the **cylindrical tower** or its extended gripper arm! This will damage the instrument!

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1. Cleaning Spills

- Turn off the instrument.
- Wipe the spill using an absorbent material. Always pull the liquid to the front of the instrument, away from the control panel and around the bath positions.
- Wipe the deck dry with a dust-free towel and 70% ethanol (or similar).
- Dispose of the material in accordance with good laboratory practices.

2. Disposing of Reagents

Chemicals used on this instrument may be hazardous. Adhere to good laboratory practices when disposing of any hazardous chemicals.

G. Compliance

1. Electromagnetic Compatibility (CE)

All instruments shipped to the European Union (EU; formerly known as the European Community) have the “CE” label on the back of the instrument, signifying that these instruments comply with the Electromagnetic Compatibility and Low Voltage Directives.

2. Restriction of Hazardous Substances (RoHS)

The RoHS directive was adopted in February 2003 by the European Union. It restricts (with exceptions) the use of six hazardous materials (including lead and mercury) in the manufacture of various types of electronic equipment. All SciGene instruments carrying the CE mark are also RoHS compliant.

3. Electrical and Electronic Equipment Waste (WEEE)

WEEE is a European Community directive that became European Law in February 2003. The crossed out wheeled bin symbol on a product or its packaging indicates that it must NOT be disposed of with standard waste at the end of its lifetime. Instead, it is your responsibility to return it to a designated collection point for electronic equipment recycling. Contact your local SciGene distributor for information on drop off locations.

4. Laboratory Equipment Standards

This symbol indicates the instrument meets Laboratory Equipment standards UL 61010-1:2004 R7.05 and CAN/CSA-C22.2 61010-1:2004 as certified by TÜV Rheinland of North America, Inc.

5. Electromagnetic Compatibility (EMC)

The **Little Dipper Processor** is a Class A digital device under FCC Title 47 Part 15B and designated as Class A electrical equipment for measurement, control, and laboratory use (EN61326).

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II. UNPACKING AND SET UP

A. Unpacking the Instrument

The Little Dipper Processor weighs about 59 lbs (26.5 kg) and is generally shipped in one carton containing the instrument along with an interior box for accessories. Caution is needed when lifting the unit to protect yourself and others from personal injury. Two people are required to lift the unit, one on each side, keeping it balanced and sharing the load.

1. Open the large carton and remove the protective foam.
2. With the assistance of another person, place fingers under the bottom of the instrument and carefully lift out and place on the bench.
3. Cut the tie wrap securing the gripper arm.

NEVER lift the unit by the robot arm. Always grip from the bottom when lifting!!

If the instrument is damaged upon arrival, please retain all shipping materials and contact SciGene or your local distributor for assistance.

B. Parts Provided

The following items may be included with the instrument or in separate carton, depending on configuration:

- Touchscreen with cable
- Power cord
- Digital thermometer
- Centrifuge buckets
- Baths
- Slide racks
- Pads, spacers or covers
- User Manual


C. Environmental Requirements

Ensure that the area where the instrument is installed meets the following conditions, for reasons of safety and performance:

Ambient temperature	+15 to +32°C (58 to 90°F)
Relative humidity	20 to 80% RH non-condensing
Air flow clearance	6 inches (15 cm) minimum at back of unit

D. Installation and Location

Place the instrument on a lab bench or stable, level surface of at least 30 x 22 inches (76 x 56 cm) with a vertical clearance of 30 inches (76 cm) within a few feet of a power source that delivers 20 Amps of constant power. Position the instrument with 6 inches of clearance at the back to allow adequate air flow of the cooling fan and ensure unimpeded access to the main power and emergency shut off (EMO) switches on the right side of the instrument.

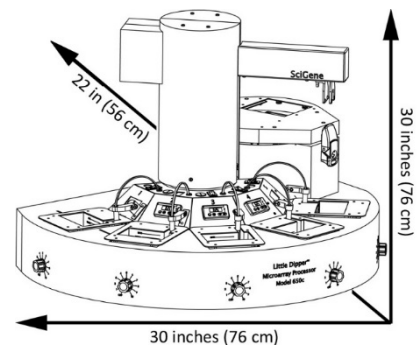E. Installing Baths

1. Clean baths, slide racks and stir bars by rinsing in running tap water, followed by 100 ml of DI water and then 100 ml of ethanol. Dry with a lint-free towel.
2. Clean temperature sensors with DI water and a lint-free towel.
3. Insert baths in the black basins. Rotate temperature sensors up prior to inserting any standard or low volume heatable (LVH) baths, and then ensure all temperature sensors are lowered down prior to starting a protocol.
4. Place a stir bar in each bath.

Never pour any liquid into the black basins! A removable bath must be inserted prior to adding buffer.

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Never heat a bath without buffer! Baths may need more buffer added due to evaporation.

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F. Installing Centrifuge Bucket Assemblies


Always use the correctly sized centrifuge buckets, spacers and pads matching your sample rack: 24-position (large) or 12 position (small).

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
To install a centrifuge bucket assembly:

1. Unlock and open the centrifuge door.
2. Gently squeeze the side tabs of the **red** bucket and slide it onto the pegs of the rotator arms marked "BALANCE RACK".
3. Gently squeeze the side tabs of the **green** bucket and slide it onto the pegs of the rotator arm marked "SAMPLE RACK".
4. Place an absorbent pad in each bucket then a black frame spacer on top.
5. Place a slide rack into the **red** balance bucket with an equal number of slides as those to be processed.
6. Close and lock the centrifuge.

24-Slide Rack Assembly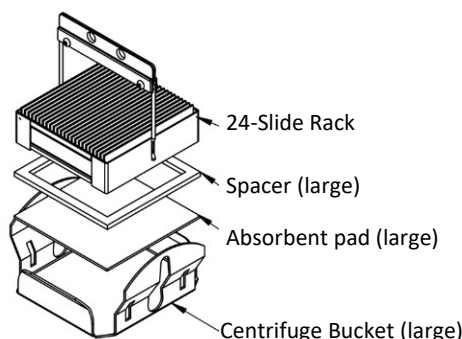


12-Slide Rack Assembly

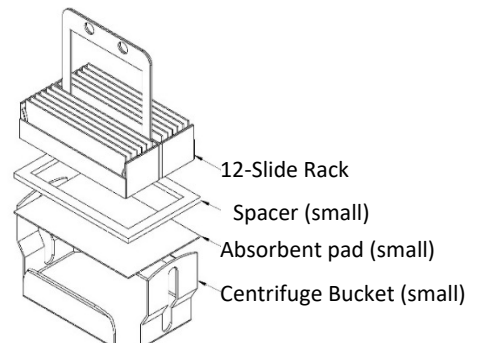III. USING YOUR LITTLE DIPPER PROCESSOR


A. Components, Controls and Accessories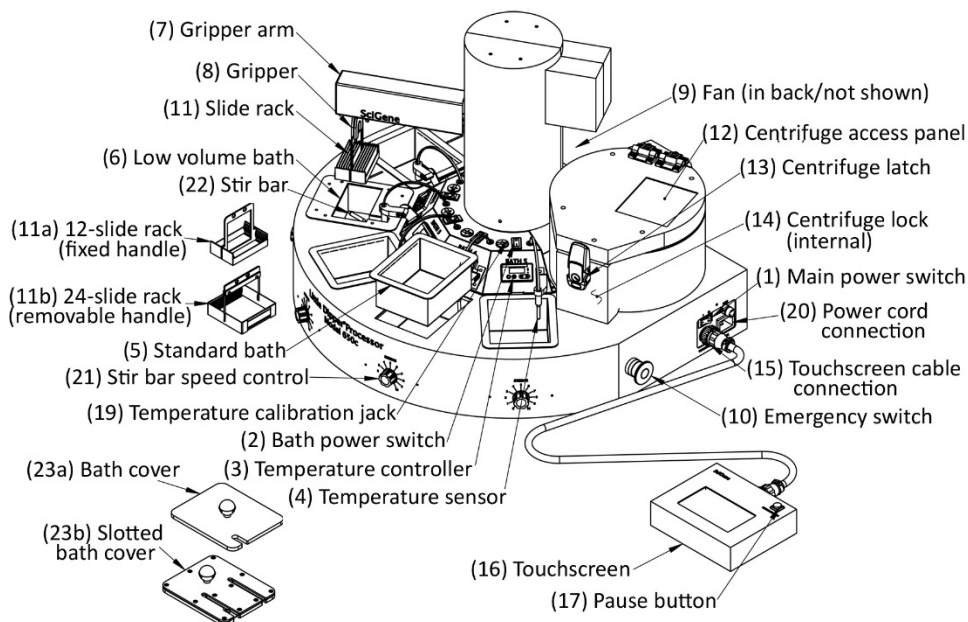

Name	Function
Main power switch (1)	Turns on power to the unit
Bath power switch (2)	Turns on power to each bath, for heating liquid
Temperature controller (3)	Used to set and observe bath temperature
Temperature sensor (4)	Transmits bath temperature to controller.
Standard bath (5)	Removable insert, compatible with most liquids/solvents
Low volume bath (6)	Removable stainless steel insert
Gripper arm (7)	Rotates and moves slide racks up and down in baths
Gripper (8)	Opens, closes and secures the slide rack assembly
Fan (9)	Draws air through the instrument cabinet for cooling
Emergency switch (10)	Cuts power for safety
Slide rack (11a,b)	Holds slides during processing.
Centrifuge panel / trap door (12)	Automatically opens for the gripper to insert a slide rack
Centrifuge latch (13)	Keeps lid closed
Centrifuge lock (14)	Internal lock for safety
Touchscreen cable connection (15)	Receptacle for attaching control cable
Touchscreen (16)	Used to create, store and run protocols and for calibration
Pause button (17)	Toggle switch that stops and restarts a running protocol
Temperature calibration jack (19)	Receptacle for cable to thermometer
Power cord connection (20)	Receptacle for power cord
Stir bar speed control (21)	Rotating switch to adjust stir bar speed
Stir bar (22)	Creates a gentle vortex in the bath
Bath cover (23a)	Lids help achieve and maintain temperatures above 70°C

**Optional accessories*


Internal View of Centrifuge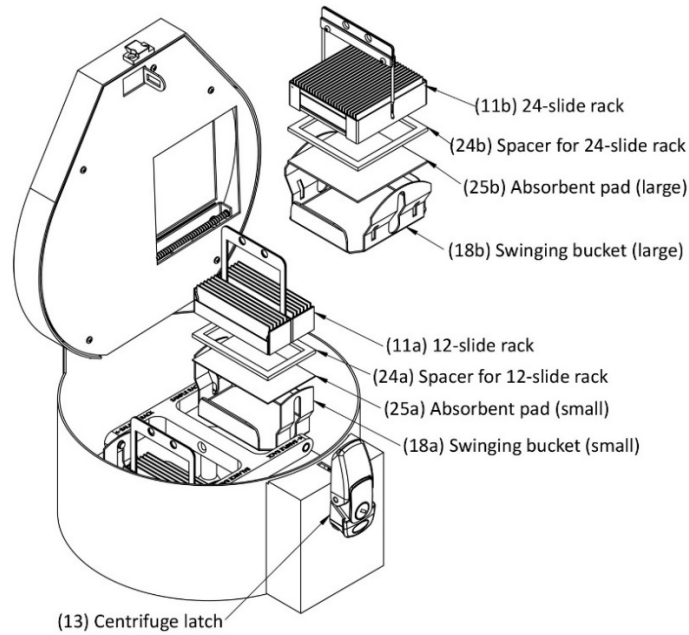

Name	Function
Slide rack (11b)	Holds slides during processing.
Centrifuge latch (13)	Keeps lid closed
Centrifuge swinging buckets (18a,b)	Green (sample) and red (balance) rack buckets
Slide rack spacer (24a,b)	Raises rack above absorbent pad
Absorbent pad (25a,b)	Captures liquid beneath slide rack during centrifuge spin


B. Touch Screen and Power Connections

1. Connect the braided cable to the touchscreen and to the receptacle labeled "TOUCHSCREEN" on the right side of the instrument, using the circular connectors.
2. Connect the power cord to the receptacle labeled "POWER" on the right side of the instrument and to a power source that delivers 20 Amps or more for 115V units or 10 Amps or more for 230V units. Avoid connecting to a circuit with any other equipment that draws significant power.
3. Turn on the main power switch on the right side of the instrument. The touchscreen will illuminate and motors will activate.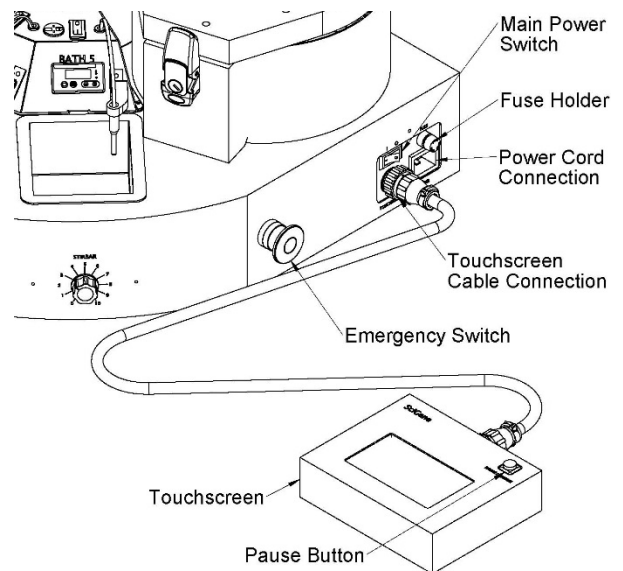

C. Using the Temperature Controller

The Little Dipper Processor is equipped with a Watlow EZ-Zone temperature controller calibrated at the factory to provide accurate temperatures from 35 to 95°C. The controller has two LED displays and four push buttons. The left display shows ACTUAL temperature in red. The right display shows the SET or programmed temperature in green. To enter the SET temperature, simply push the up and down arrows until the desired temperature is shown in the green display on the right. The instrument will adjust the heat until the SET temperature is attained.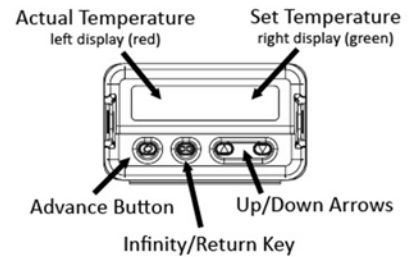

D. Calibrating the Temperature Controller

The EZ-Zone controllers are calibrated at the factory to provide accurate bath temperatures from ambient 35 to 95°C. Re-calibration is recommended *only if* the display varies by more than one degree (1°C) to the temperature shown on an externally connected NIST certified digital thermometer (Cat. #1051-52-0).

To adjust a controller to achieve accurate temperatures:

1. Turn ON the instrument and set the controller to 65°C. Allow 30 minutes for temperature to stabilize.
2. Using the cable provided with the digital thermometer, plug one end into the blue receptacle found above the controller.
3. Turn on the thermometer. Allow 1 minute for thermometer to stabilize. The actual temperature of the bath will be displayed.
4. Calculate the difference between the temperature shown on the controller and on the digital thermometer. For example, if the controller displays 65.0°C and the thermometer displays 66.5°C, the difference (offset) is 1.5°C.
5. On the controller, press the up and down arrows simultaneously for 3 seconds. The left display shows "A1" and the right display shows "open".
6. Press the Advance Button (green circle) 3 times until the right display shows "i.CA". The left display shows the offset value between the controller and thermometer when the unit was last calibrated.
7. Use the up or down arrow to adjust the offset value to the temperature difference calculated in Step 4. For example, if the controller shows a temperature 1.5°C lower than the thermometer, adjust the offset by adding 1.5 to the value shown on the display.
8. Press the Infinity Key (∞) twice to exit calibration and return to the operation display. Verify that the temperature on the thermometer matches the display. The controller is now calibrated to provide accurate temperatures from 35 to 95°C.


IV. SYSTEM OPERATION

A. Using the Touchscreen

Use the handheld touchscreen to navigate between screens, select and run protocols, create new protocols and control instrument movement.


B. Using the Pause Button

Instrument movement can be stopped and resumed during a protocol using the red “PAUSE/RESUME” button at top right on the touchscreen. When pressed, the instrument will stop and the button light will flash. To resume the protocol, press pause again.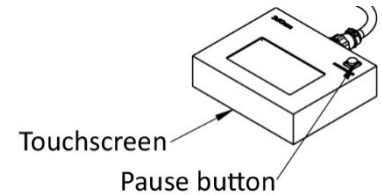

C. Using the Emergency Stop (EMO)

To immediately stop movement of the gripper arm and stir bars, press the red Emergency Stop (EMO) button on the right side of the instrument. Bath heaters will continue to receive power but automated movement will cease and touch screen will turn off. **Restore power to the gripper arm and stir bars by rotating the EMO button a quarter turn clockwise.**

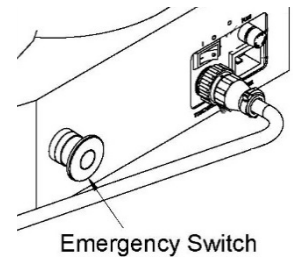If the Emergency off switch is used to stop the instrument, please allow 30 seconds for the touchscreen to properly shut down before turning it on.

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Press to stop a protocol.

Twist clockwise to resume.

The protocol will NOT resume upon releasing the EMO switch and must be restarted.

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D. Using a Bath Cover for Pre-Heating

A bath cover (SciGene cat. #1080-12-0) may be used to speed pre-heating (up to 95°C) and reduce buffer evaporation, prior to slide processing.

Slide a cover over the bath to achieve high temperatures and reduce evaporation.

Bath covers must be removed before starting any protocol unless the protocol includes a User Controlled Pause in the step prior to the bath with the protocol.

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E. Loading a Slide Rack

The pre-loaded **QUIK TST** protocol provides the operator an opportunity to practice loading the slide rack onto the gripper. An empty slide rack is moved automatically from bath to bath with up/down agitation, followed by brief centrifugation. No liquid is required. Power switches for baths should remain in the OFF position.

1. Ensure that baths and centrifuge buckets are properly installed and correspond to size of the slide rack to be used.
2. Place an empty slide rack in the **red** balance bucket in the centrifuge.
3. Verify all temperature sensors are rotated down and power to all baths are turned off.
4. Turn on main power located on the right side of the instrument.
5. Using the touchscreen, press **Run Protocol** from the home screen and select **QUIK TST**.
6. Hold an empty slide rack in your left hand centered in Bath #1 and aligned with the black lines on the instrument as shown below.
7. Press **Run Protocol** on the touchscreen. The arm will lower into Bath #1 with the gripper open. When the gripper arm reaches the end of its downward travel, align the holes in the handle with the two bumps in the gripper paddle as it slowly closes.


Hold slide rack in Bath #1 to load.

Place and hold rack on gripper bumps.

Release rack as gripper closes.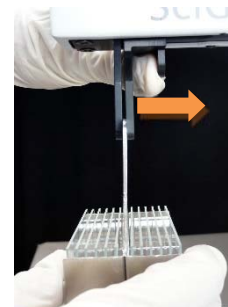

Use thumbpad to manually open gripper.

If the slide rack does NOT load properly, press the pause button and use the thumbpads on the gripper to manually open and load the rack. Press pause again to continue.

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8. The instrument will agitate the rack in each bath and then transfer it to the centrifuge. When the protocol is complete, unlock and open the centrifuge lid to remove the rack.

F. Running a Programmed Protocol

The instrument comes configured with validated protocols installed. Refer to the Little Dipper Processor Method corresponding to your protocol for details on how to set up the instrument. Method sheets can be found online at www.scigene.com.

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The following are general guidelines on how to set up and run a pre-installed protocol.

1. Ensure baths, racks and stir bars are clean and free of buffer residue and dust.
2. Insert baths and stir bars as specified for the protocol.
3. Rotate all temperature sensors down.
4. Add recommended buffer(s) to the fill line on the side of the baths.
5. Turn on main power to the instrument.
6. Turn on power to bath(s) to be heated and set temperature(s). Use a standard bath cover for temperatures above 70°C to reduce heating time and evaporation. Allow time for baths to reach set temperature(s).
7. Open the lid of the centrifuge and place a slide rack with the same number of slides to be processed into the **red** balance bucket.
8. Add the slides to be processed into the sample rack.
9. Remove any bath covers.
10. Submerge the slide rack in Bath #1 using your left hand between the two black lines.
11. On the touchscreen home page, press **Run Protocol**, then select the desired protocol and press **Run Protocol**.
12. When the gripper arm stops in the down position in bath #1, align the holes of the slide rack handle onto the bumps gripper paddles and wait for the gripper to close. The instrument will now perform the steps as programmed.
13. At completion, remove the rack from the centrifuge. Some protocols end with the rack remaining in the gripper. Open the gripper to release the slide rack by pushing the thumb tab on the rear gripper paddle.
14. Clean the instrument per section **VI. C. Cleaning After Every Use**.

V. USING CUSTOM PROTOCOLS


A. Creating or Editing a Protocol

Custom protocols can be created using the touchscreen. All protocols start from Bath #1.


1. From the touchscreen Home page, press Create/Edit Protocol. On the **Create/Edit Protocol** screen, touch <NEW> on the screen.

If editing a protocol, select it directly from the Protocols list.


2. Begin by highlighting Step 1 of the protocol. On Step 1, press the green button under each parameter to input the value for the step. Descriptions of the available parameters are below.
3. Press Save Step after finishing the inputs for each individual step. If you move to another step without saving, that line will be lost.
4. Following the input of all the steps, press Save/Save As. This will pull up a keyboard to enter the protocol name.
5. Press Exit to return to the Home page.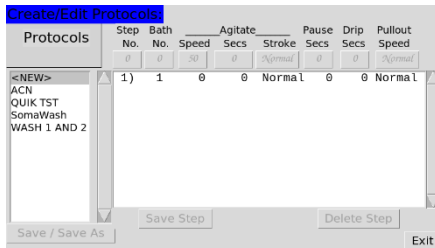


Highlight <NEW> on the Create/Edit Protocols page to begin creating a new protocol.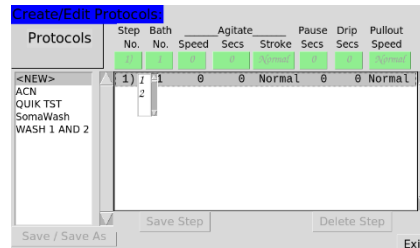

Highlight line 1 to begin input of the parameters and use the pull down menus to enter the Step No., Bath No., Agitate Speed, Agitate Stroke, and Pullout Speed.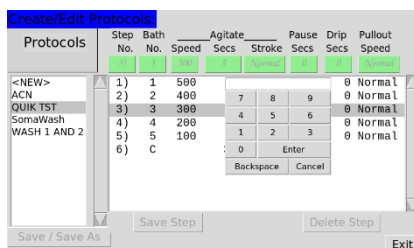

Use the touchpad to enter the number of seconds for the Agitate, Pause, and Drip functions.

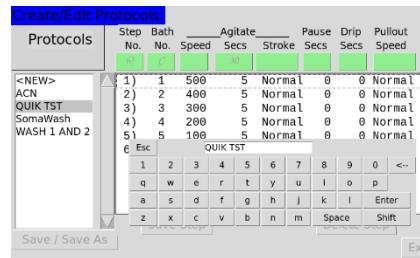Press Save/Save As and use the keyboard to name the protocol. Press Enter to Save or Esc to go back and continue editing.

Available Parameters

Step No.:	Choose from the pull-down menu.
Bath No.:	Choose from positions 1-5 or Centrifuge from the pull-down menu. Step 1 must always be Bath 1. The Centrifuge may only be the final step.
Agitation Speed:	Choose 0-600 from the pull-down menu.
Agitation Secs:	Input the number of seconds from the touchpad for the step.
Agitation Stroke:	From the pull-down menu, choose <i>Normal</i> or <i>Long</i> . A <i>Normal</i> stroke agitates below the surface of the liquid. A <i>Long</i> stroke brings the slides above the surface of the liquid each stroke.
Pause Secs:	Input the time in seconds for the rack to remain in the bath following agitation before moving to the next step.
Drip Secs:	Enter a time to hold the rack above the bath before the next step.
Pullout Speed:	Choose between a <i>Normal</i> and <i>Slow</i> withdraw from the bath. A slow withdraw is recommended for the final step if the protocol doesn't utilize the centrifuge.

B. Deleting a Protocol

1. To delete a protocol, press **Control Panel** from the home screen then **Advanced Settings**.
2. Press the **Delete Protocol** button to open the screen.
3. Choose a protocol to delete then press "Delete Protocol?" Choose OK to confirm.
4. Press the BACK button on each of the previous screens to return to the home screen.

VI. MAINTAINING THE LITTLE DIPPER PROCESSOR

A. Powering Off

Turn the power switch to the OFF position and unplug the power cord before performing any service procedure. Once the instrument has been powered down, allow 30 seconds before restarting.

B. Cleaning Before Every Use

Baths, racks and stir bars must be free of buffer residue and dust before processing slides. If they were not cleaned and stored properly after the last use, wash and dry them as follows:

1. Fill baths containing the rack(s) and stir bars with tap water.
2. Wipe racks, stir bars and baths with a lint-free towel.
3. Rinse in running tap water.
4. Rinse with approximately 100-200 ml of DI water.
5. Rinse with 100 ml of ethanol and dry with a lint-free towel.
6. Clean temperature sensors with DI water and a lint-free towel.

C. Cleaning After Every Use

1. Wash baths, stir bars and racks with running tap water then rinse in DI water and ethanol and dry with a lint-free towel. Store in clean plastic bags to protect from dust.
2. Clean temperature sensors with DI water and lint-free towel.
3. Wipe exterior of instrument cabinet with a damp paper towel.

D. Periodic Cleaning and Calibration

1. Calibrate bath temperatures using a NIST-certified thermometer. See section **VI. A. Temperature Controller Calibration**.
2. Remove and wash the centrifuge buckets and spacers using warm water and lint-free towels.
3. Replace the absorbent pad in each centrifuge bucket.
4. Clean all exterior surfaces of the cabinet, gripper paddles and gripper arm housing with damp, lint-free towels.
5. Remove the centrifuge buckets. Clean the inside surfaces of the centrifuge and rotor with damp, lint-free towels.
6. Inspect the centrifuge buckets for signs of wear or deformation where they hang from the rotor. Replace if worn or bent.

VII. TROUBLESHOOTING

Symptom	Cause	Solution
Instrument does not turn on.	Blown fuse(s)	Replace fuse(s) near main power switch.
Touch screen does not illuminate when main power is on.	Not connected or EMO is engaged. / Touchscreen did not fully shutdown.	Check connections and rotate EMO switch to release the emergency stop. Cycle the power on then off . Leave off for 30 seconds then restart.
Touch screen does not respond when buttons are pressed.	Pause button is engaged or touch screen is faulty.	Press to release Pause button or contact SciGene or your local distributor for assistance.

VIII. SPECIFICATIONS

Electrical	
Voltage	115V AC; 50/60 Hz; 20A or 230V AC; 50/60 Hz; 10A
Dimensions	
Size (HxWxD)	20 x 28 x 22 inches (51x 71 x 56 cm)
Weight	59 lbs (26.5 kg) net
Performance and Controls	
Temperature Range	Ambient +5°C to 95°C
Temperature Regulation	± 0.5°C from set point
Agitation Speed	0 to 600 relative units
Centrifuge RPM/Force	690 RPM / 50 g
Digital Thermometer Output	T-type Thermocouple

IX. ORDERING INFORMATION

Cat. #	Description	UoM
1080-40-1	Little Dipper® Processor for Agilent, 115V	EA
1080-40-2	Little Dipper® Processor for Agilent, 230V	EA
1080-55-1	Little Dipper® Processor for Somalogic, 115V	EA
1080-55-2	Little Dipper® Processor for Somalogic, 230V	EA
1080-70-1	Little Dipper® Processor for FISH, 115V	EA
1080-70-2	Little Dipper® Processor for FISH, 230V	EA
1080-10-1	Standard Bath, 670 ml.	EA
1080-10-5	Low Volume Heatable Bath, 275 ml.	EA
1080-12-0	Bath cover for Little Dipper. Use for pre-heating only.	EA
1080-20-1	Slide Rack, 12-Position for standard 3 inch slides. Includes handle.	EA
1080-20-5	Slide Rack, 24-Position for standard 3 inch slides. Includes handle.	EA
1080-21-0	Absorbent Pads for Slide Rack Buckets, 4-Position and 24-Position, 25/PK.	PK
1080-21-1	Absorbent Pads for Slide Rack Buckets, 12-Position, 25/PK.	PK
2022-00-2	CytoZyme Stabilized Pepsin, 50X Concentrate, 20 ml. Store at 4°C.	EA
2022-00-3	CytoZyme Stabilized Pepsin, 50X Concentrate, 100 ml. Store at 4°C.	EA
2022-10-2	CytoZyme Reaction Buffer, 1L. Store at RT.	EA
2022-10-3	CytoZyme Reaction Buffer, 4L. Store at RT.	EA
2010-00-1	FISH Wash Buffer 1 (0.4xSSC/0.3% IGEPAL, pH 7), 4L. Store at RT.	EA
2010-00-2	FISH Wash Buffer 2 (2xSSC/0.1% IGEPAL, pH 7), 4L. Store at RT.	EA

X. DECLARATION OF CONFORMITY

Little Dipper® Processor

SciGene
 3350 Scott Blvd., Bldg. 39-02
 Santa Clara, CA 95054 USA

Declares that the above referenced product(s) meets the essential requirements of the following European Union Directives by using the relevant standards shown below to indicate compliance.

EMC Directive 2004/108/EC

Standards:

EN 61326-1	2006	Electrical equipment for measurement, control and laboratory use to include:
EN 61000-4-2	1995	1998 +A2: 2001
EN 61000-4-3	2006	
EN 61000-4-4	2004	
EN 61000-4-6	2007	

LVD Directive 2006/95/EC

EN 61010-1	2010	Safety requirements for measurement, control and laboratory use 3rd Edition Part 1: General requirements
------------	------	---

IEC 60204-1	2006	Electrical equipment of machines
-------------	------	----------------------------------

RoHS Directive 2011/65/EU

Terry Gill

 Name of Authorized Representative

Sunnyvale, California, USA

 Place of Issue

Director of Product Manufacturing

 Title of Authorized Representative

February 13, 2023

 Date of Issue

 Signature of Authorized Representative